


AquaClean

Microbiocidal Additive for Water Soluble Heating Bath Fluids



Water Baths

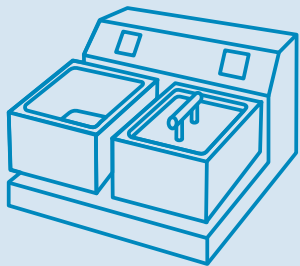

AquaClean inhibits growth of bacteria, algae, fungi and precipitation of inorganic salts.

AquaClean shows a blue indicator color. Bacterial contamination is indicated by decoloration of the fluid.

AquaClean does not contain aldehydes or azides.

Application:

- Add 5ml AquaClean to 1 liter of heating bath fluid.
- Renew fluid and additive every 4 weeks. In case the blue color of fluid has faded, renew earlier.

Compatible for use as an additive in water reservoirs inside cell culture incubators. (Its non-volatile active ingredients cannot invade cell-cultures via air.)

Cat.#	WAK-AQ-250-50L 250 ml concentrate sufficient for 50 liters of fluid	Active ingredients:	Surfactants, quaternary ammonium compounds, complexing agents
Precautions:	Avoid eyes and skin contact	Storage:	Room temperature
Disposal:	Biologically degradable	Materials compatibility:	Universal except for aluminium and acrylic glass

scientifix

PO Box 662 South Yarra
VIC 3141 Australia

www.scientifix.com.au

Free call: 1800 007 900
T: +61(0)3 8540 5900
F: +61(0)3 9543 7827
E: info@scientifix.com.au

**WAK - Chemie
Medical GmbH**

